


GrainSense

Technical sheet

The handheld GrainSense device measures the quality of cereal grains and other crops in seconds: **protein, moisture, carbohydrates, and oil contents***

* Percentages are calculated on a dry, wet or fixed basis, based on country guidelines.

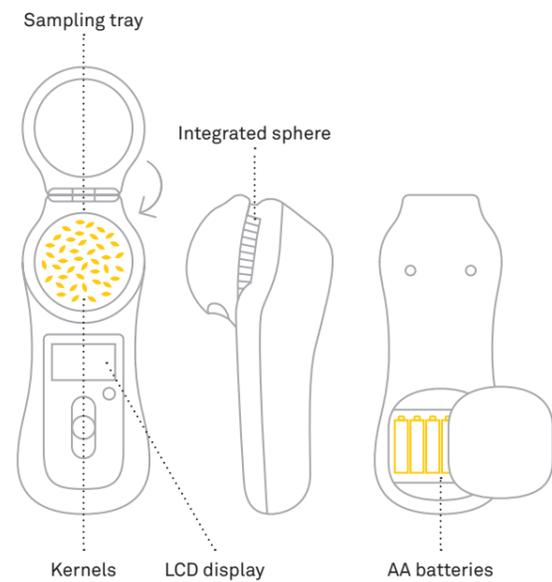

The technical principle is **Near-infrared (NIR) spectroscopy** in the so-called third overtone wavelength range. This technique has been used in laboratory instruments for years. GrainSense is the first to realize such an instrument in a handheld format.

Because of the **patented sampling technology** (grain inside an “integrating sphere”) the light intensity arriving at the detector is several hundred times higher than otherwise possible. This enables the building of a **small, battery-operated device**.




GrainSense
360° light penetration method (integration sphere)

- + Works with simpler and more affordable technology
- + Device can be small
- + Enables small samples
- + Wider use than grain as other types of samples possible
- + Short measurements time

Technical specifications

| | |
|------------------------|---|
| Size | Hand-held (footprint 270 mm x 115 mm) |
| Weight | 820 grams (without batteries) |
| Batteries | 6 x AA Alkaline (recommended to use batteries designed for industrial/heavy use) |
| Battery operation | 50 to 150 measurements depending on battery quality and type of use |
| Measurement principle | Near infrared transmittance spectroscopy |
| Sample size | ≈ 3 grams (60-80 cereal kernels) |
| Measurement time | About 30 seconds, including the device warm-up and the user loading the sample |
| Species | E.g. wheat, barley, rye, oat and rapeseed New species can be added according to needs / Note: Species are country specific |
| Operational conditions | +5 to +45 C 20 to 90 % RH (non condensing) |
| Storage temperature | -10 to +60 C |
| Protection | Designed for outdoor use - except raindrops on the sample tray will affect the moisture result |
| Bluetooth | LE 4.1 |
| Language | Multiple languages (based on ISO 8859-1 character set) |
| Mobile application | Android/iOS |

The key components of the GrainSense solution are the GrainSense device, mobile application, and cloud-based database:

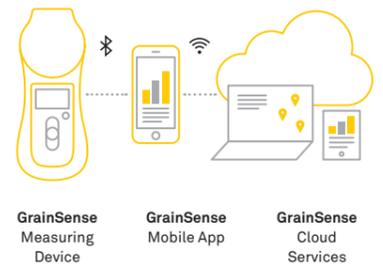

1. GrainSense Device: measures and analyzes the sample and interacts with the GrainSense Mobile application via Bluetooth.

2. Mobile application: connects with cloud account and downloads calibrations and other settings to the Device and uploads measurements results to the database.

3. Cloud based storage: to backup the measurement results.

Customer support: support@grainsense.com

